

HA0325T - TFA2 and PSY240 – LED Lighting Strip - Fitting Instructions

This information sheet details instructions for fitting the LED Lighting Strip.

For the continuing safety of the operators of the equipment and for the protection of the equipment itself, the operator should take note of the **Warnings, Cautions and Notes** throughout this leaflet, and, where applicable, in any associated literature.

1.1 Installation

- 1) Remove all packaging.

Item (2) is applicable only if fitting to the Nexus Table Frame Overhead Shelf unit.

- 2) Using the M6 hex bolts supplied, attach the spring clips to the underside of the shelf frame.

Item (3) is applicable only if fitting to something other than the Nexus Table Frame Overhead Shelf unit.

- 3) Discard the M6 hex bolts supplied, and using other suitable screws/bolts attach the spring clips to the surface where the lighting strip is to be installed.

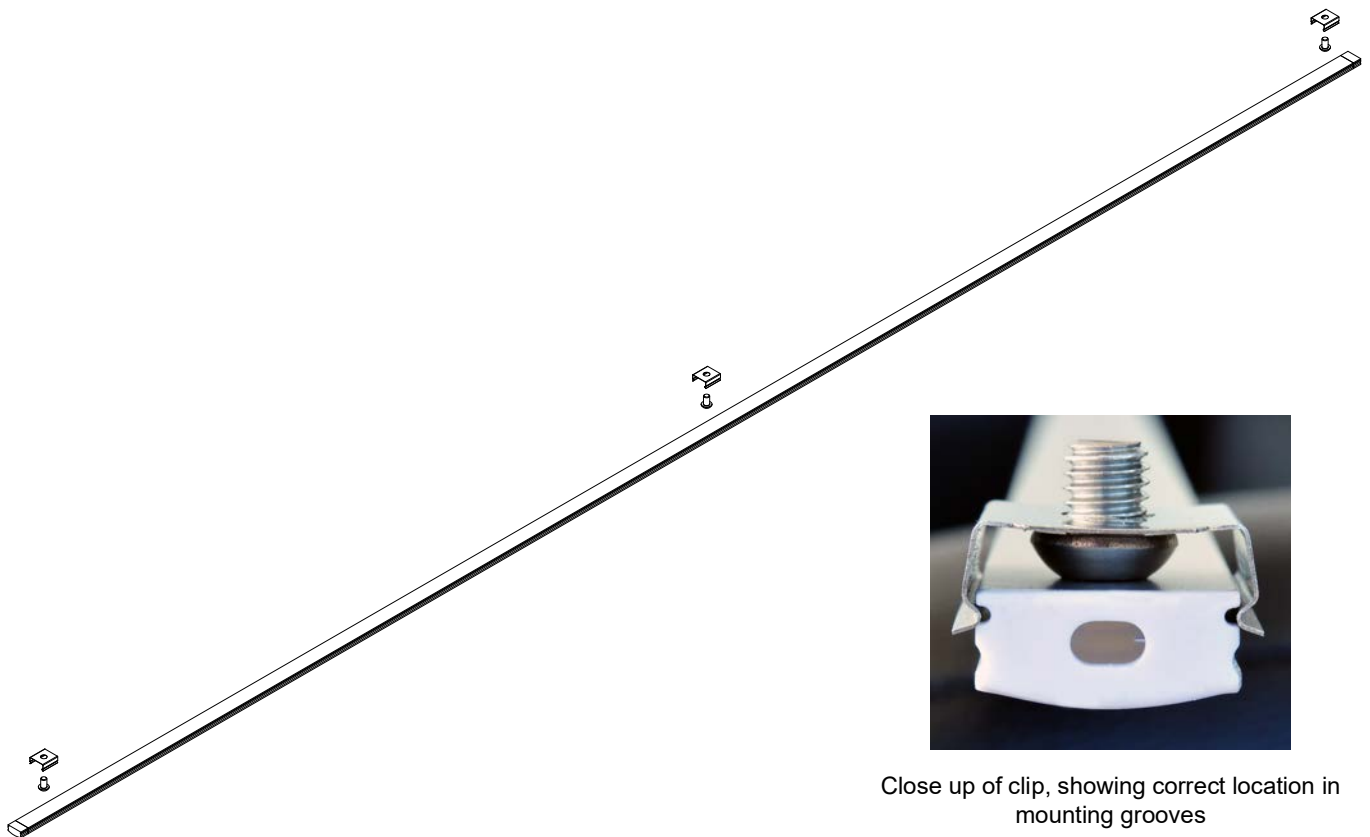


Fig. 1.1 Installing the LED light strip

- 4) Fit the lighting strip into the spring clips, ensuring that the clips locate correctly into the groove along the edges of the strip as shown above.

Thorlabs Worldwide Contacts

For technical support or sales inquiries, please visit us at www.thorlabs.com/contact for our most up-to-date contact information.



USA, Canada, and South America

Thorlabs, Inc.
sales@thorlabs.com
techsupport@thorlabs.com

Europe

Thorlabs GmbH
europe@thorlabs.com

France

Thorlabs SAS
sales.fr@thorlabs.com

Japan

Thorlabs Japan Inc.
sales@thorlabs.jp

UK and Ireland

Thorlabs Ltd.
sales@uk.thorlabs.com
techsupport.uk@thorlabs.com

Scandinavia

Thorlabs Sweden AB
scandinavia@thorlabs.com

Brazil

Thorlabs Vendas de Fotônicos Ltda.
brasil@thorlabs.com

China

Thorlabs China
chinasales@thorlabs.com



THORLABS
www.thorlabs.com